

# ***Life Diagnostics, Inc.***

610-431-7707 • Fax: 610-431-7818

Info@lifediagnostics.com • www.lifediagnostics.com

## **PEG Blocking Buffer**

Catalog No.	PEGB125-1
Volume	125 ml
Composition	Proprietary PEG-free buffer containing BSA as blocker and a proprietary disaccharide as stabilizer
Instructions	We recommend that 96-well plates be coated overnight using 100 $\mu$ l/well of antigen containing bicarbonate buffer. Wash with distilled or de-ionized water and block with 150 $\mu$ l/well PEGB125-1. Incubate for 30 minutes at room temp, remove blocker and either vacuum or air-dry. Plates should be stored desiccated.
Storage	2-8°C

Comments: Blocking and stabilizing buffer for use in ELISA or western-blot detection of PEG, PEGylated proteins and anti-PEG antibodies. This product can be used together with our PEG-free dilution buffer (PEGD50-1) and wash buffer (PEGW50-20).

Rev 062013

P.O. Box 5205 • West Chester PA 19380 • USA